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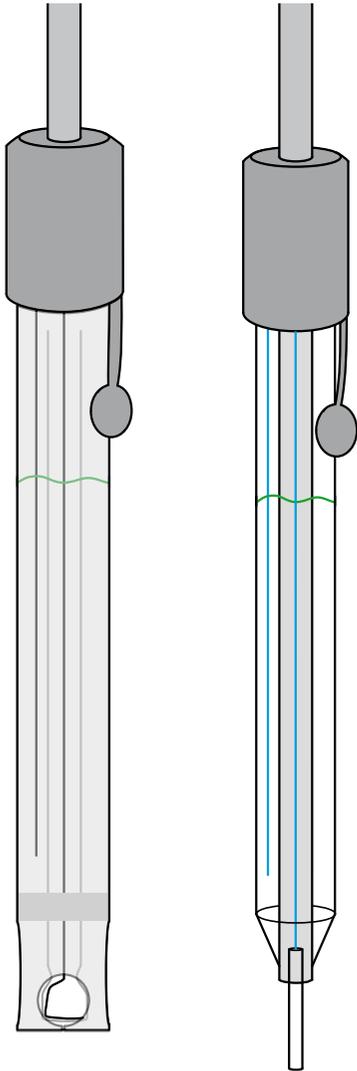
Thank you for purchasing an **ORP combination electrode or argentometry combination electrode**.

For the best results with your electrode:

- read this user manual carefully,
- comply with the precautions for use.

The ORP electrode is used to measure a solution's oxidation reduction potential.
ORP = Oxidation Reduction Potential.

The argentometry electrode is used to measure the concentration of silver ions in a solution. These electrodes may contain mercury.

 These products are highly poisonous, even at low doses. They may have very varied physical effects: nausea, vomiting, headache, loss of consciousness or other more significant problems leading to death.

 These products may be included in one or more of the following categories: carcinogenic, mutagenic, toxic for human reproduction. They may also modify the functioning of certain organs (liver, nervous system), attack the lungs and cause allergies (asthma).

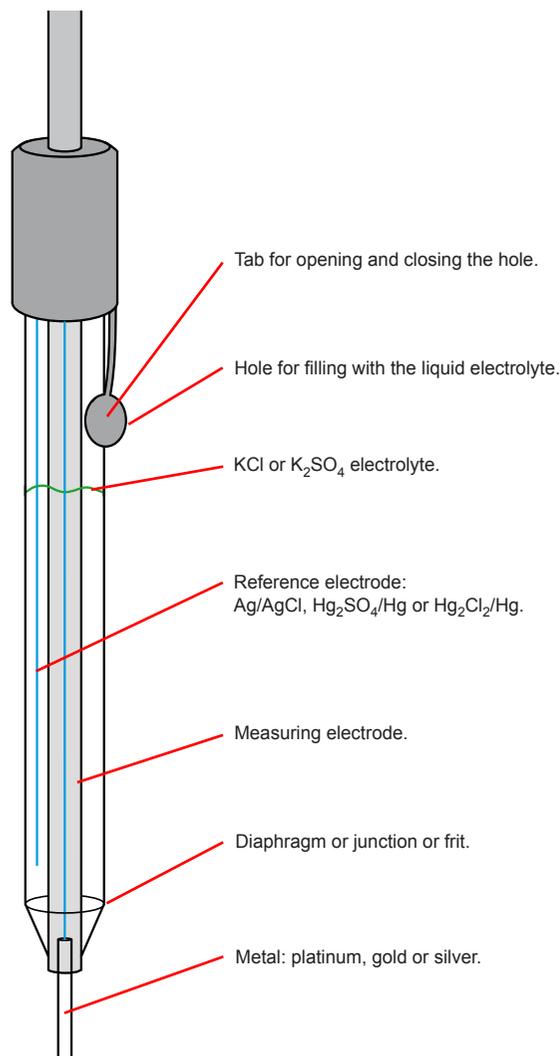
 These products may have harmful effects on the environment, in particular on organisms in aquatic environments: fish, crustaceans, algae and other aquatic plants.

 The CE marking indicates compliance with the European directives, particularly LV and EMC.

 The crossed-out dustbin means that, in the European Union, the product is the subject of selective disposal as per directive DEEE 2002/96/EC: this equipment must not be treated as household waste.

USE

- Remove the storage reservoir cap.
- If you notice the presence of an air bubble at the level of the glass membrane or diaphragm, remove it by shaking the electrode.
- Connect the electrode to the measuring instrument.
- Open the opening tab.
- Calibrate the electrode with a reference standard solution and then rinse it.
- Immerse the tip of the electrode in the solution to be measured, while keeping it vertical and making sure that the metal and diaphragm are immersed, but without exceeding the level of the internal electrolyte.
- Wait for the measurement to stabilize.
- Rinse the electrode with demineralized water after each measurement.
- Close the tab again when you have finished measuring.

The presence of white crystalline deposits on the electrode is a normal phenomenon due to evaporation of the storage solution in the reservoir cap or the internal electrolyte. To remove this deposit, simply rinse the electrode with demineralized water and then wipe it off with a clean, lint-free cloth.

Do not use aggressive and/or abrasive products (hydrofluoric acid, acetone, etc.).

Contamination of the diaphragm is the most frequent cause of measurement problems.

- Remove any impurities deposited on the glass surface of the electrode and on the outside of the diaphragm with demineralized water or with a soft cloth if the deposit is significant
- If rinsing is not sufficient, it is possible to use various chemical cleaning solutions, depending on the type of contamination:

Type of clogging	Electrodes concerned	Cleaning solution
By silver sulphide	Electrode filled with a KCl electrolyte	Solution of thiourea in hydrochloric acid
By silver chloride		Concentrated ammoniac solution
By proteins	Any electrode with an electrolytic compartment	Solution of pepsin + HCl
Other types of clogging		Ultrasonic bath with water or HCl 0.1 mol/L

MAINTENANCE

Keep your electrode in a vertical position in a reservoir cap containing an electrolytic solution identical to the electrode's internal electrolyte (KCl 1 mol/L in most of our electrodes). The level of the storage electrolyte must be lower than that of the internal electrolyte with which the electrode is filled.

Electrode reference	Filling electrolyte	Storage solution
BRPt1	KCl 1 mol/L	KCl 1 mol/L storage solution
XRPt1		
BRAG1	Saturated K ₂ SO ₄	K ₂ SO ₄ storage solution

For electrodes with liquid electrolytes, close the tab of the filling hole before storing.

Regularly check the condition of your electrode, the cleanness of the diaphragm and metal and the level of the electrolyte.

If the level of the electrolyte in the electrode seems insufficient or contaminated, you will have to top it up or empty it and refill it with a new electrolyte solution suitable for your electrode.

STATE AT DELIVERY

The ORP combination electrode or argentometry combination electrode is delivered in a cardboard box with:

- a shim,
- a storage reservoir cap,
- a bilingual user's manual,
- a multilingual accompanying datasheet.

For accessories and replacement parts, please visit our website:

www.chauvin-arnoux.com

LIFE SPAN AND WARRANTY

ORP electrodes and argentometry electrodes are wear parts. The life span of your electrode depends on the conditions in which it is used and the maintenance which you carry out on it.

The warranty does not apply if the glass is broken.

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